*vacuubrand*

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Instructions for repair: 999211-02 / version 10.01.2011  
Valid for Diaphragm pumps NT and NT VARIO  
Vacuum gauges CVC 3000 and DCP 3000

It is the users' responsibility to check the validity of the pump specific drawing and spare parts list on the basis of the serial number of the pump.  
If the pump is older or newer please contact VACUUBRAND and order the pump specific drawing and spare parts list stating the pump type, serial number and year of manufacture.

Kundendienst und Service:  
After sales service: Telefon: 09342/808-209  
Contact your local dealer or call (++49) 9342/808-193.

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## **! DANGER**

- Danger! Immediate danger. Death or severe injuries as well as damage to equipment and environment can occur.

## **! WARNING**

- ☞ Warning! Possible danger. Severe injuries as well as damage to equipment and environment can occur.

## **! CAUTION**

- Caution! Possible danger. Slight injuries as well as damage to equipment and environment can occur.

## **NOTICE**

Note. Disregarding of notes may cause damage to the product.

Caution! Hot surface!

Isolate equipment from mains before removing the cover.

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## Safety information!

### NOTICE

Ensure that repair is done only by an appropriately trained and supervised technician. Obey local and national safety requirements. Ensure that the technician is familiar with the safety procedures which relate to the products processed by the pumping system. Ensure that the pump is decontaminated before repair and that you take adequate precautions to protect people from the effects of hazardous substances if contamination has occurred.

Read the **instructions for use** for the equipment carefully, especially sections "Safety informations" and "Use and operation". If necessary order "Instructions for use" from VACUUBRAND.

### CAUTION

**Wear parts have to be replaced regularly.** In case of normal wear the lifetime of the diaphragms and valves is > 10000 operating hours. Bearings have a typical durability of 40000 h. Motor capacitors have a typical durability in the range of 10000 to 40000 h depending strongly on the operation conditions like ambient temperature, humidity or load.

- Check every capacitor regularly by measuring its capacity and estimating its operation time. Exchange old capacitors early enough to prevent a failure. If an overaged motor capacitor fails it might get hot and even melt and may cause a flame to form which could be **dangerous for persons and equipment in the vicinity**. The capacitors have to be replaced by an electrician.

### NOTICE

If hazardous or corrosive substances have been pumped:

- The device will be contaminated with the process chemicals which have been pumps during operation. Ensure that the pump is decontaminated before maintenance.
- Take adequate precautions to protect people from the effects of dangerous substances. Wear appropriate safety-clothing, do not inhale and avoid contact with skin.

Before starting repair check that the required parts are available and of the correct type before you start your work.

- Check the operating sequence mentally using exploded view drawings, spare parts lists and circuit diagram on feasibility, safety requirements and consequences on safety and function of the equipment.

### WARNING

**- Isolate equipment from mains** and **wait two minutes** before starting maintenance to allow the capacitors to discharge.

**- Vent the pump, isolate the pump and other components from the vacuum system.** Allow sufficient cooling of the pump. Drain condensate, if applicable.

**☞ Ensure that the pump cannot be operated accidentally. Never operate the pump if covers or other parts of the pump are disassembled. Never operate a defective or damaged pump.**

Ensure that **maintenance** is done only by suitably trained and supervised technicians. Ensure that the maintenance technician is familiar with the safety procedures which relate to the products processed by the pumping system.

### CAUTION

In order to comply with law (occupational, health and safety regulations, safety at work law and regulations for environmental protection) vacuum pumps, components and measuring instruments returned to the manufacturer can be repaired only when certain procedures (see section "**Notes on return to the factory**") are followed.

## **DANGER**

- Use only **genuine spare parts and accessories**. Otherwise safety and performance of the equipment as well as the electromagnetic compatibility of the equipment might be reduced.  
Possibly the CE mark or the C/US conformity becomes void if not using genuine spare parts.

## **CAUTION**

- Replace damaged parts in any case.

Check operability and safety after repair.

Check the electrical safety (protective conductor resistance, insulating resistance, high voltage test) according to IEC 61010 and national regulations.

## **NOTICE**

## **CAUTION**

- In case of problems return the pump to the factory for inspection or repair if necessary.

Obey regulations when disposing of solvents and chemicals.

In order to comply with law (occupational, health and safety regulations, safety at work law and regulations for environmental protection) vacuum pumps, components and measuring instruments returned to the manufacturer can be repaired only when certain procedures (see section "Notes on return to the factory") are followed.

### Scraping and waste disposal

Dispose of the pump and any components removed from it safely in accordance with all local and national safety and environmental requirements.

Particular care must be taken with components and waste oil which have been contaminated with dangerous substances from the process.

- You may authorize us to dispose the equipment at your expense.

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## General notes

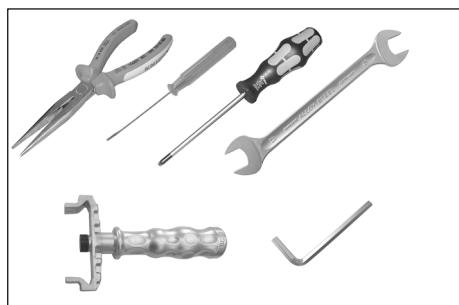
Inspection and maintenance of the pump heads and **replacement of diaphragms and valves**: See instructions for use of the diaphragm pump.

All parts of the pump head can be replaced individually.

### Note regarding the set of disc springs and undular spring (636277)

Depending on type and version diaphragm pumps may contain a set of disc springs or a special undular spring to prestress the bearing on the shaft. Replace defective parts only by parts of the same type.

**Tools** required (details for all types of pumps):

- Phillips screw driver size 2
- Open-ended wrench w/f 17 / 20
- Hex key size 5
- Diaphragm key w/f 66 (order no.: 636554)
- Torx screw driver TX20
- Screw driver 2.5 mm
- Flat pliers
- 2 big screw driver

## Replacing the fan cover and the fan wheel

Unscrew three fixing screws at the fan cover.

Remove pump feet at the lower side of the fan cover.

Remove fan cover.

Use two big Philips screw driver to lift off the fan wheel.

Remove tolerance ring from the motor shaft and assemble a new one.

Press in fan wheel on the motor shaft as far as it will go.

Position fan cover and align.

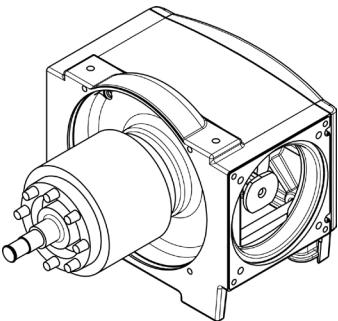
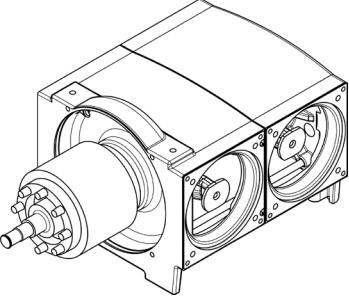
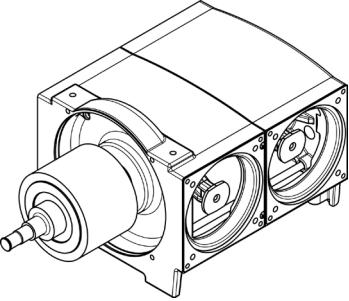
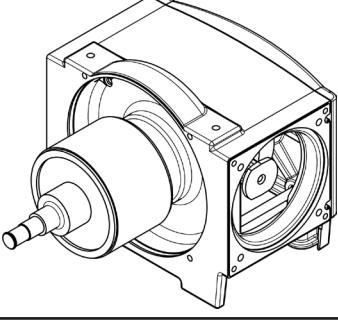
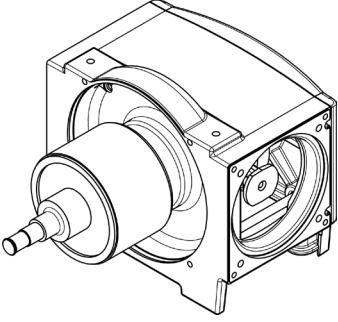
Apply a drop of adhesive (e. g. Omnifit 50M) on the thread of the fan cover fixing screws. Screw in screws. Max. torque 1.5 Nm.

Apply a drop of adhesive (e. g. Omnifit 50M) on the thread of the pump feet. Turn in pump feet.

## Replacing the pump aggregate

### NOTICE

We recommend to replace always the complete pump aggregate!

Order no. pump aggregate	suitable for	
636343	MZ 2C NT VARIO PC 3002 NT VARIO	
636342	MD 4 NT VARIO MD 4C NT VARIO MV 2 NT VARIO PC 3003 NT VARIO PC 3004 NT VARIO	
636344	ME 8 NT ME 8C NT MD 4NT MD 4C NT MV 2 NT PC 600 series	
636345	MZ 2D NT	
636346	MZ 2 NT MZ 2C NT ME 4 NT ME 4C NT ME 4R NT PC 500 series	

Disassemble the pump from the pumping unit if necessary.

Disassemble pump heads (see instructions for use, section „Replacing diaphragms and valves“).

Disassemble handle.

Unscrew three fixing screws at the fan cover.

Remove pump feet at the lower side of the fan cover.

Remove fan cover.

Use two big Philips screw driver to lift off the fan wheel.

Unscrew four motor fixing screws and remove motor.

Pull off rivets which mark the orientation of the head cover (if applicable) from the defective pump aggregate and insert rivets into the corresponding positions of the new aggregate.

If the orientation of the head cover is given by a mark (notch) at the housing, apply a mark at the corresponding position of the new housing.

Position motor on the new aggregate.

Apply a drop of adhesive (e. g. Omnifit 50M) on the thread of the motor fixing screws. Screw in screws, torque 6 Nm.

Press in fan wheel on the motor shaft as far as it will go.

Position fan cover and align.

Apply a drop of adhesive (e. g. Omnifit 50M) on the thread of the fan cover fixing screws. Screw in screws. Max. torque 1.5 Nm.

Apply a drop of adhesive (e. g. Omnifit 50M) on the thread of the pump feet. Turn in pump feet.

Assemble handle.

Assemble pump heads (see instructions for use, section „Replacing diaphragms and valves“).

Assemble pump into the pumping unit if necessary.

## NOTICE

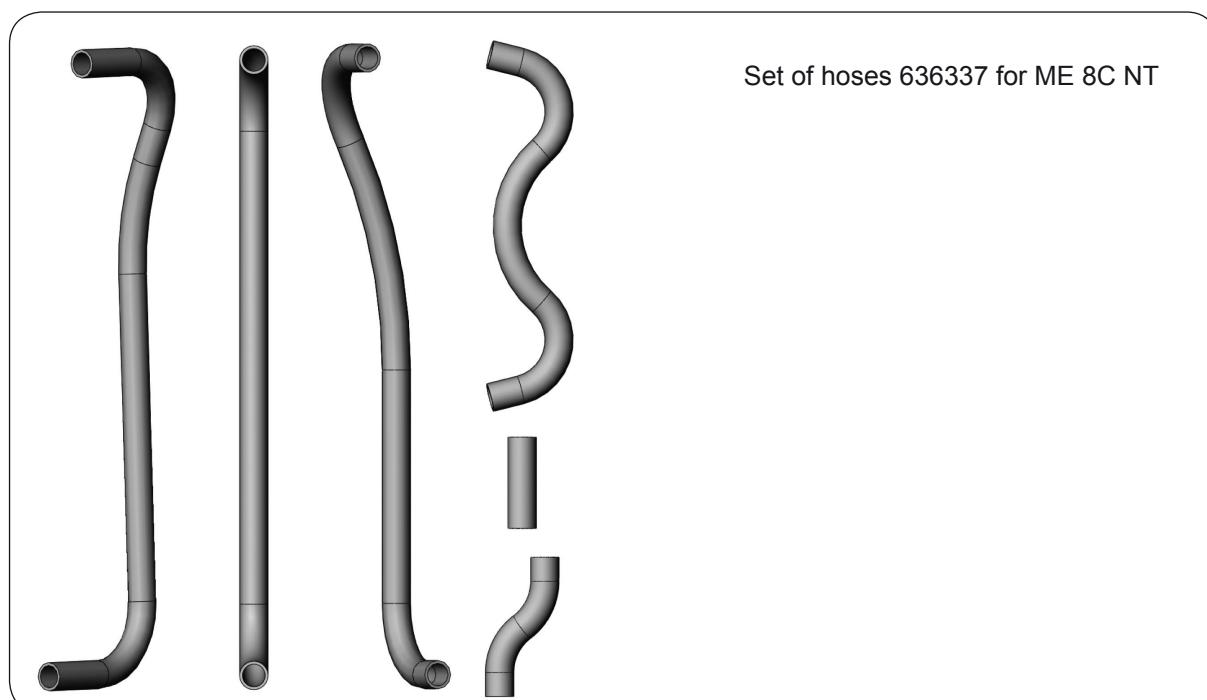
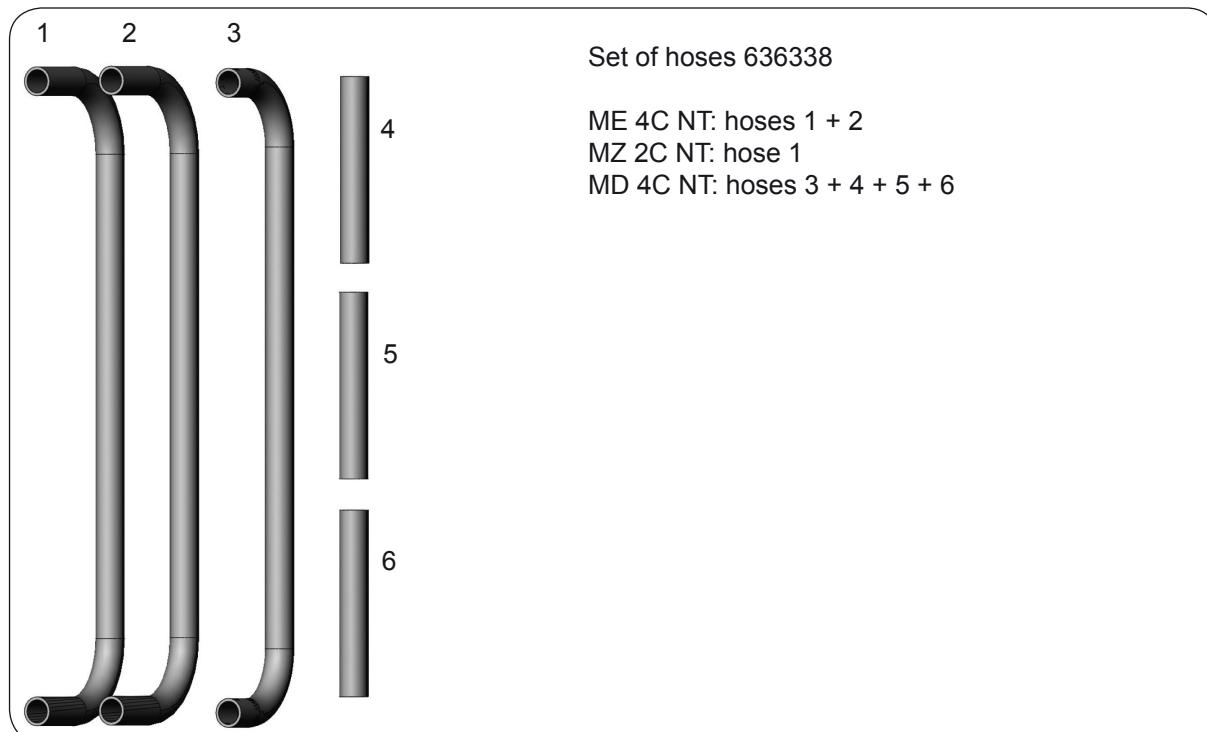
Special tools and instructions for replacing the eccentric unit with rods are available on request.

## Notes on using the sets of hoses

**Attention:** The sets of hoses are partially designated for several types of pumps and therefore the sets may contain redundant parts.

**VACUUBRAND doesn't take back redundant parts!**

Stretch the ends of the hoses before assembling by warming and using an appropriate mandrel if necessary (PTFE hoses max. 280°C).

## CVC 3000 / DCP 3000

### Replacing the front part with LCD and the pcb

Separate device from mains.

Pull off selection knob.

Unscrew the four Torx screws at the rear side of the device.

Disassemble the two housing parts carefully.

Disassemble ribbon cable from the pcb.

Unscrew the two nuts beneath the interface.

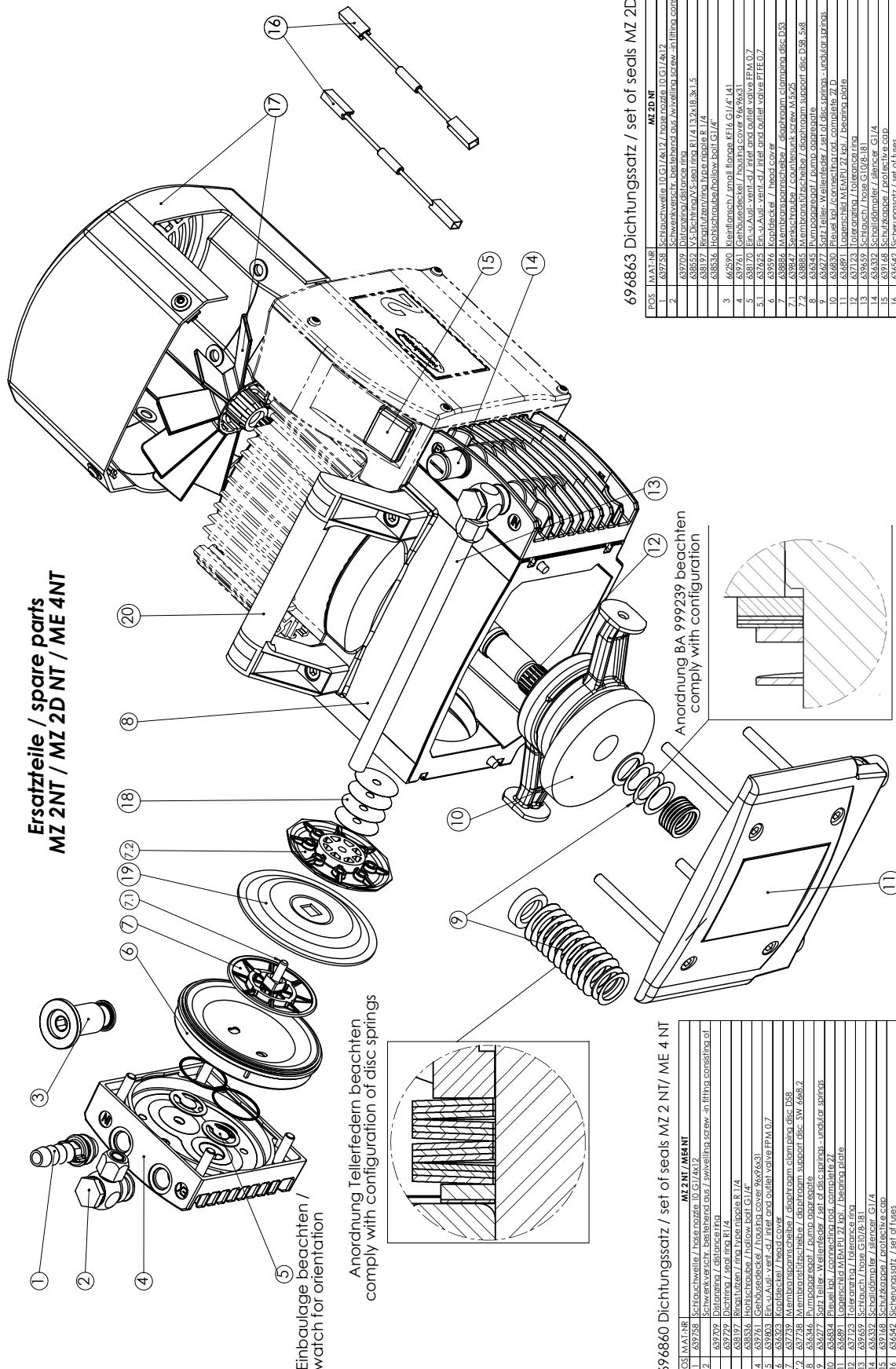
Position the new pcb in the housing rear part.

Screw in the nuts beneath the interface.


Assemble the ribbon cable of the new front part with LCD on the pcb. Comply with correct position of the ribbon cable.


Assemble the two housing parts and screw in Torx screws.

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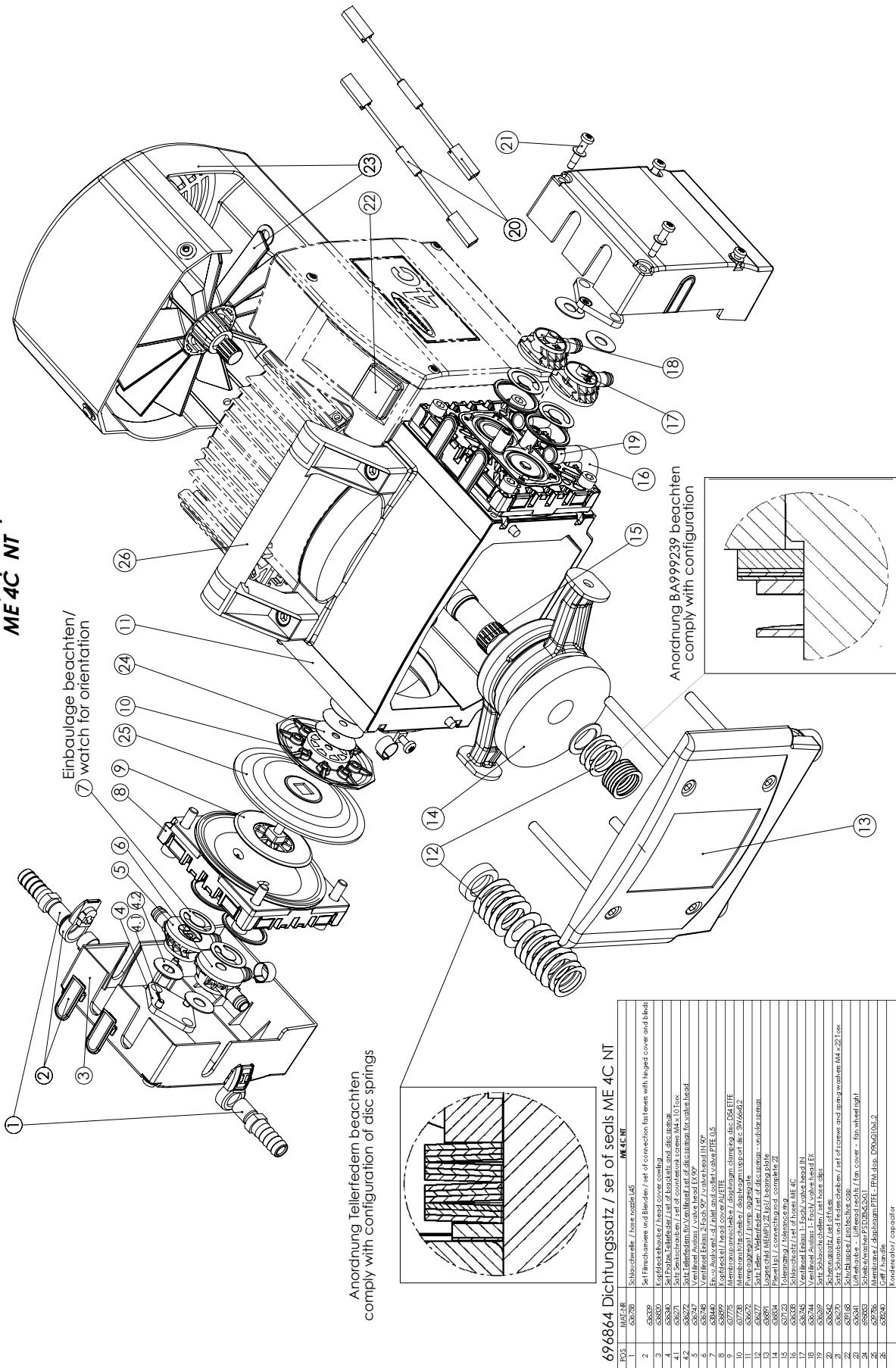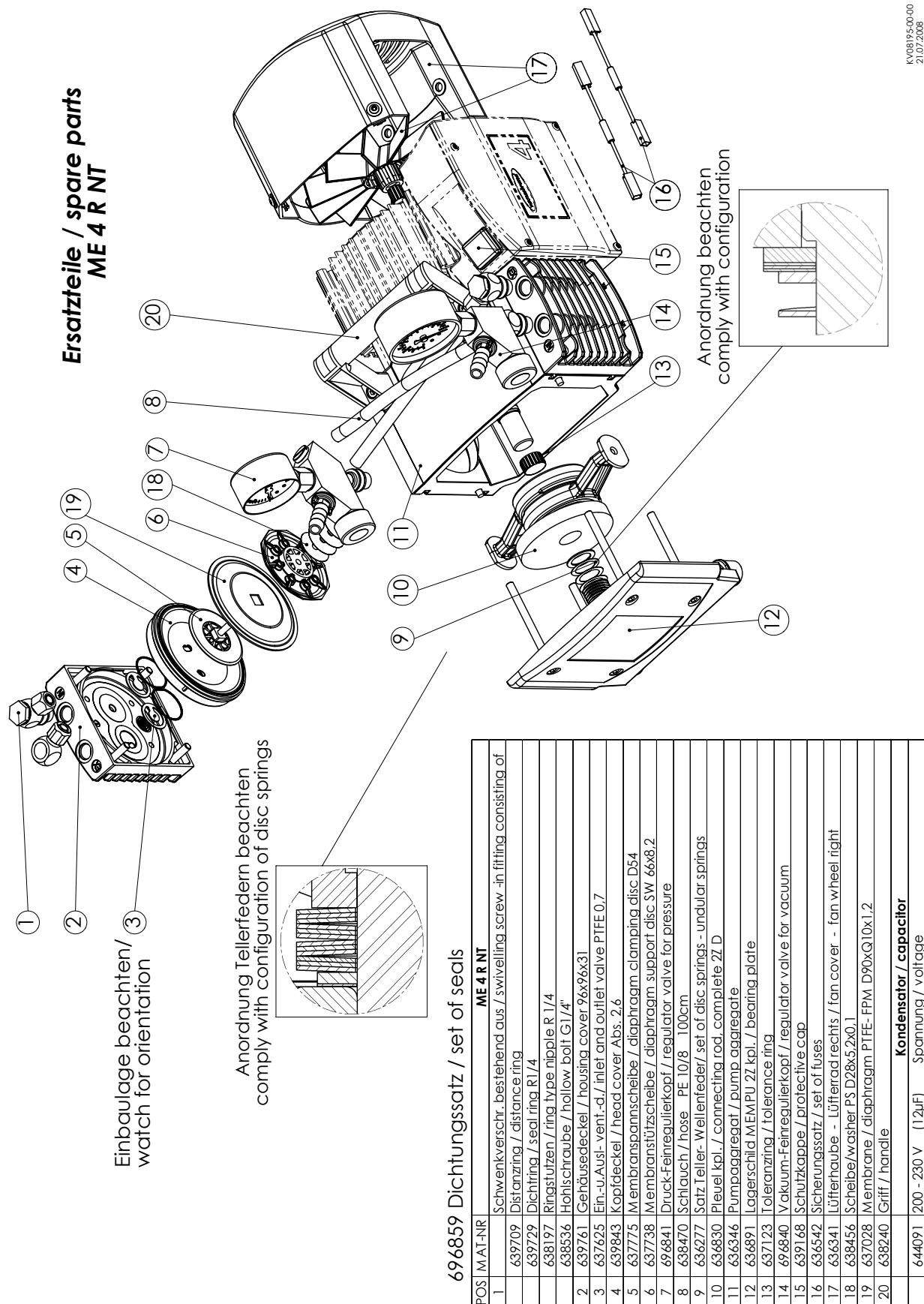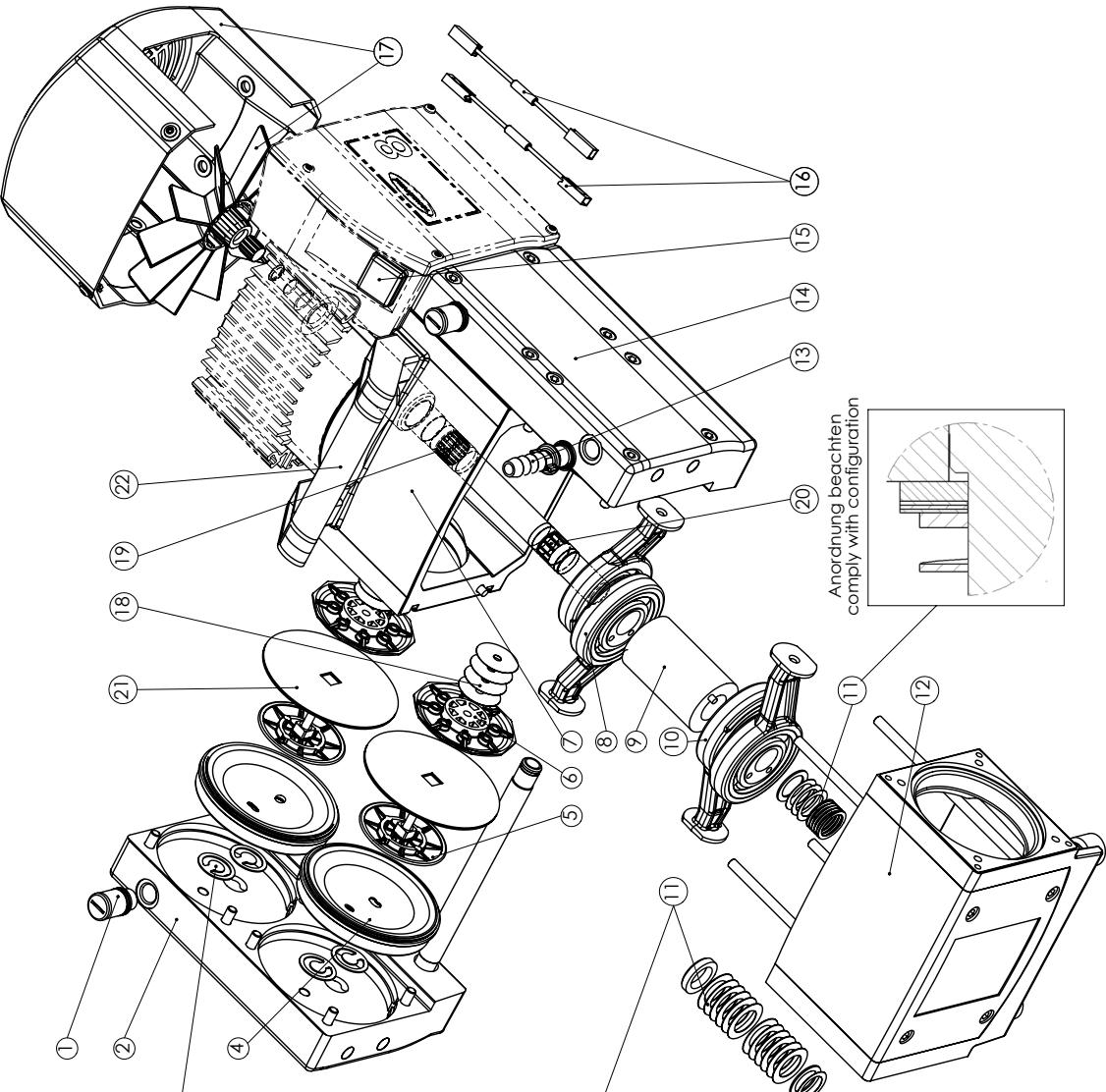
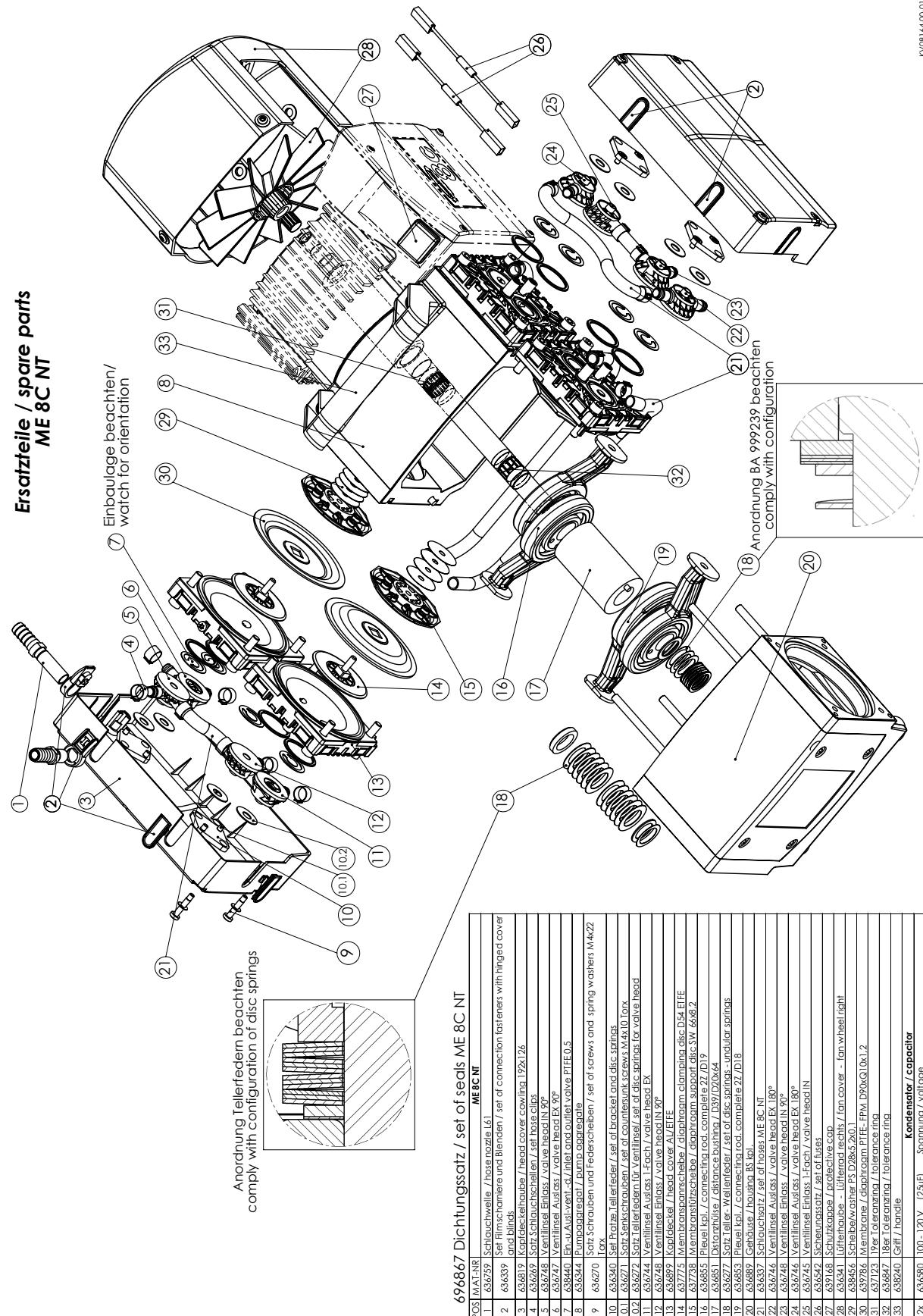
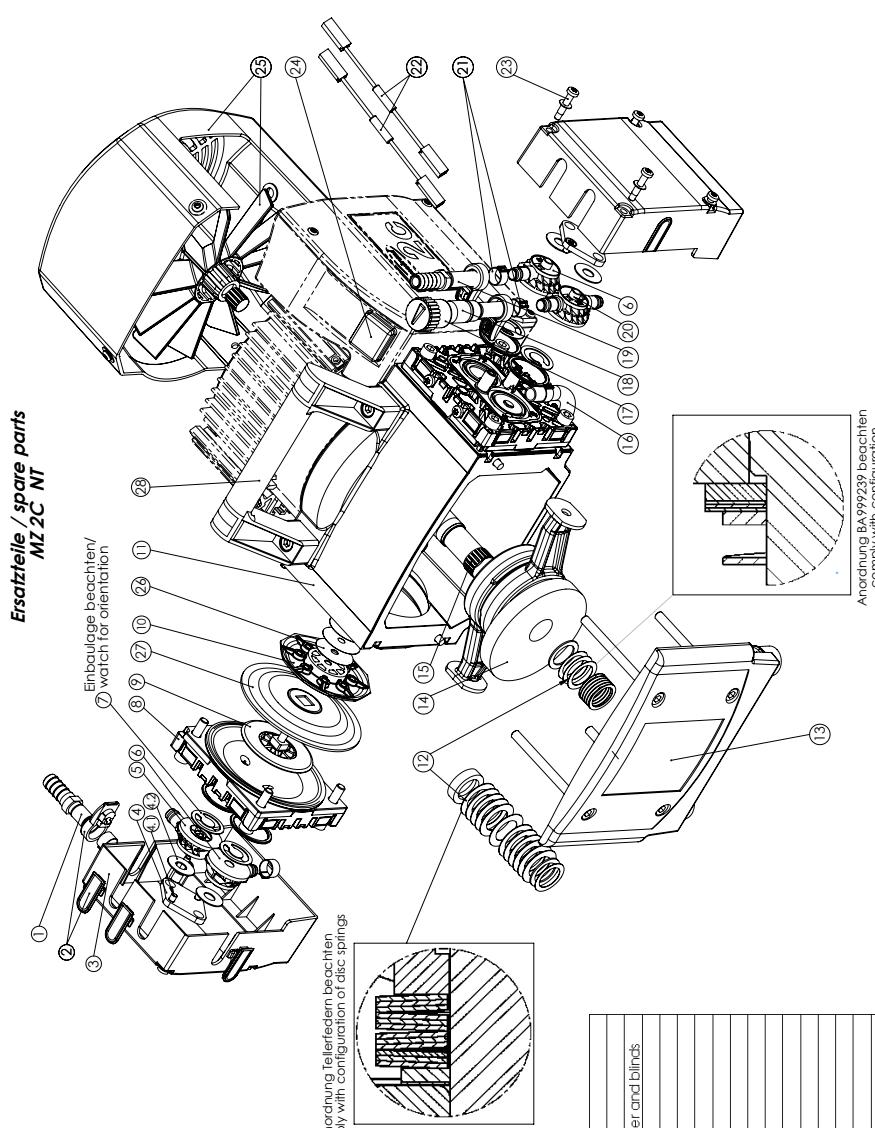
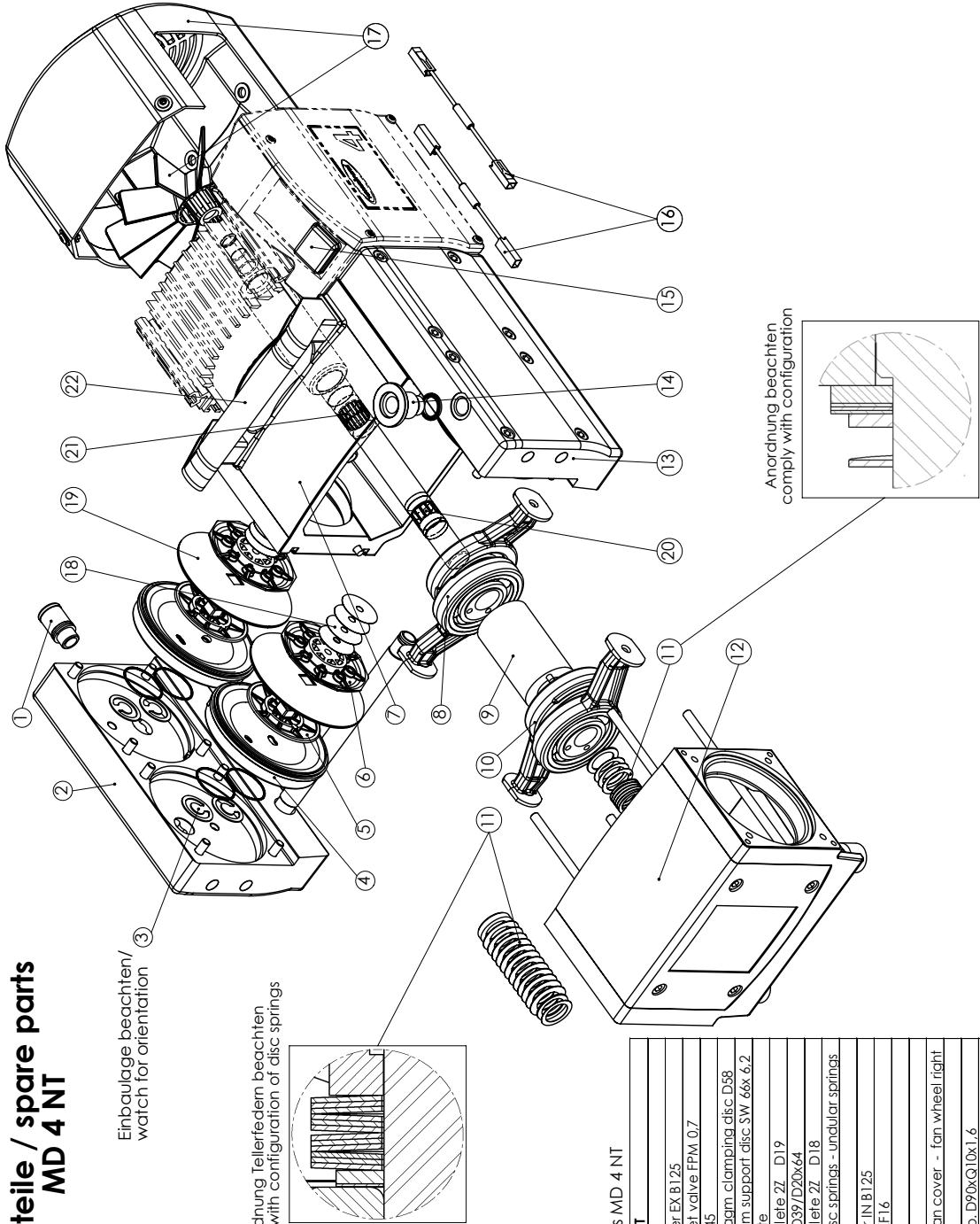
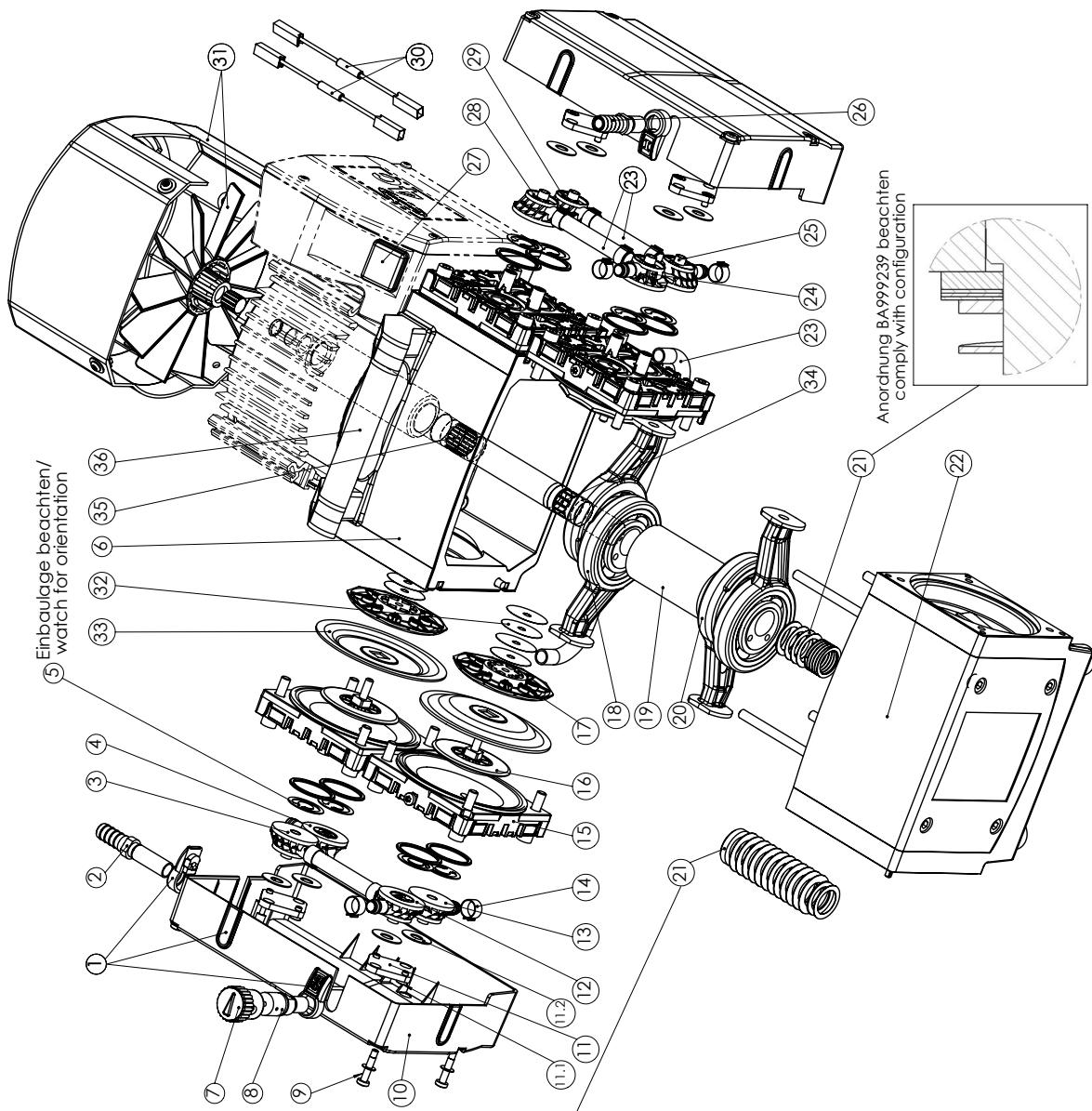
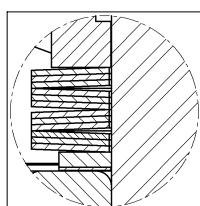
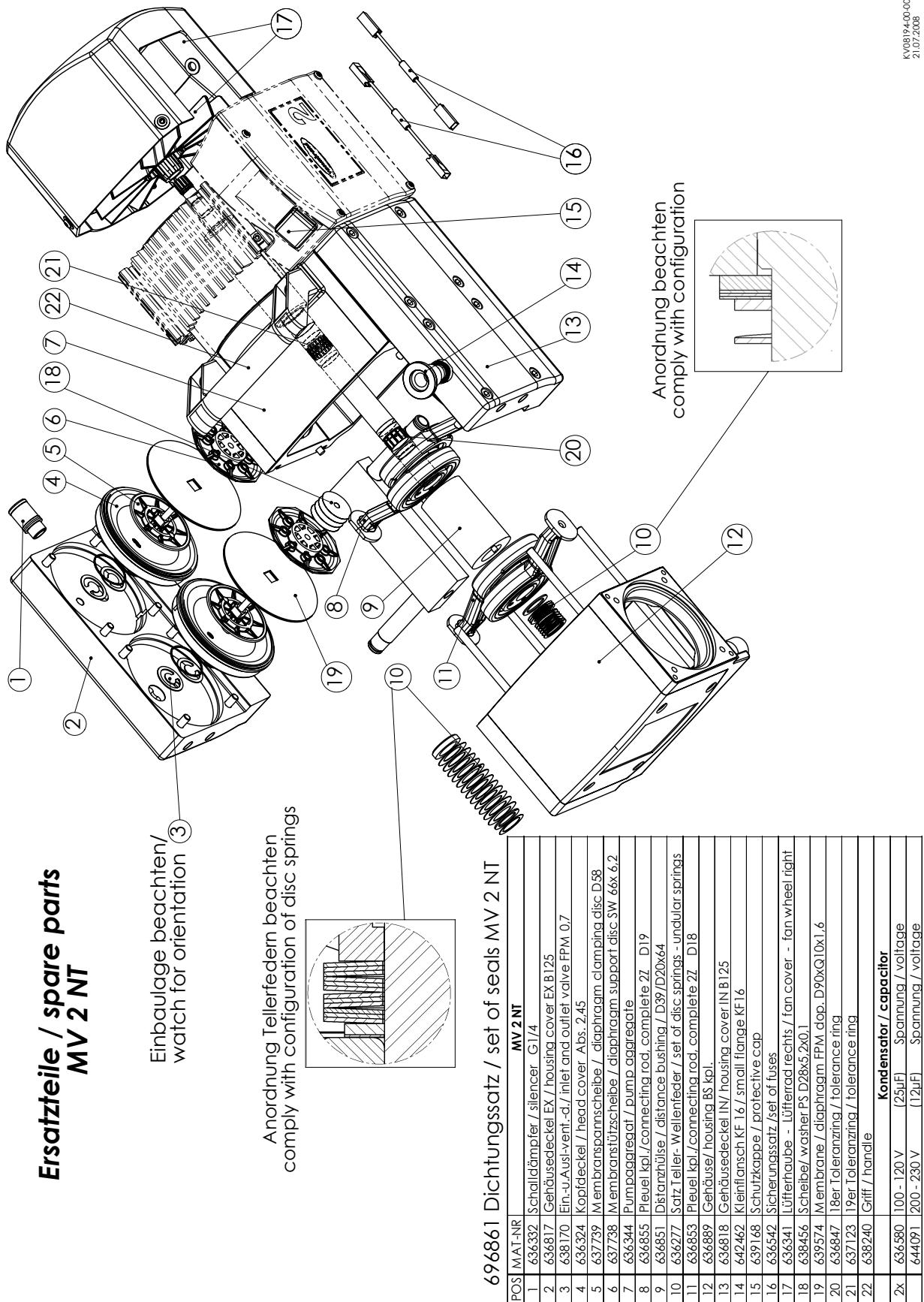
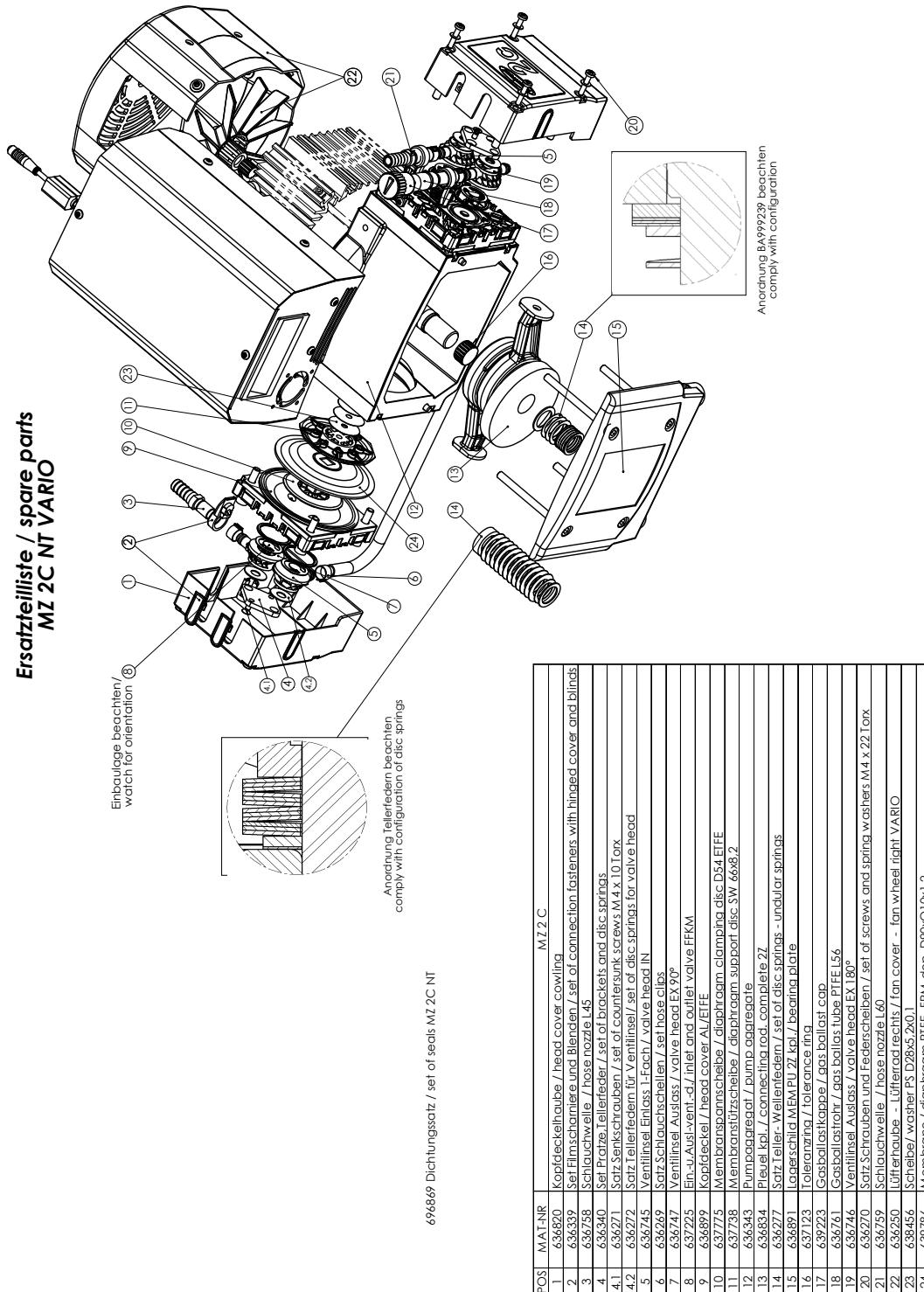
**VACUUBRAND GMBH + CO KG**

-Vakuumtechnik im System-

© 2007 VACUUBRAND GMBH + CO KG Printed in Germany

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*Documents are only to be used and distributed completely and unchanged. It is strictly the users' responsibility to check carefully the validity of this document with respect to his product. 999211 // 10/01/2011*

## Notes on return to the factory

### Repair - return - DKD calibration

#### NOTICE

Safety and health of our staff, laws and regulations regarding the handling of dangerous goods, occupational health and safety regulations and regulations regarding safe disposal of waste require that for all pumps and other products the "**Health and safety clearance form**" must be send to our office duly completed and signed before any equipment is dispatched to our premises.

Fax or post a completed copy of the health and safety clearance form to us in advance. The declaration must arrive before the equipment. Enclose a second completed copy with the product. If the equipment is contaminated you must notify the carrier.

**No repair / DKD calibration is possible unless the correctly completed form is returned. Inevitably, there will be a delay in processing the equipment if information is missing or if this procedure is not obeyed.**

If the product has come in contact with chemicals, radioactive substances or other substances dangerous to health or environment, the product must be decontaminated **prior to sending it back to the factory**.

- Return the product to us **disassembled and cleaned** and accompanied by a certificate verifying decontamination or
- Contact an industrial cleaning and **decontamination service** directly or
- Authorize us to send the product to an industrial cleaning facility **at your expense**.

To expedite repair and to reduce costs, please enclose a detailed description of the problem and the product's operating conditions with every product returned for repair. We submit **quotations** only on request and always at the customer's expense. If an order is given, the costs incurred are offset from the costs for repair or from the purchase price, if the customer prefers to buy a new product instead of repairing the defective one.

- **If you do not wish a repair on the basis of our quotation, the equipment might be returned to you disassembled and at your charge!**

In many cases, the **components must be cleaned in the factory** prior to repair. For cleaning we use an environmentally responsible water based process. Unfortunately the combined attack of elevated temperature, cleaning agent, ultrasonic treatment and mechanical stress (from pressurised water) may result in damage to the paint. Please mark in the health and safety clearance form if you wish a **repaint at your expense** just in case such a damage should occur.

We also replace parts due to optical aspects upon your request.

#### NOTICE

**Before returning** the equipment ensure that (if applicable):

- Equipment has been cleaned and/or decontaminated.
- All inlet and outlet ports have been sealed.
- Equipment has been properly packed, if necessary, please order an original packaging (costs will be charged), marked as appropriate and the carrier has been notified.
- Ensure that the completed health and safety declaration is enclosed.

We hope for your understanding for these measures, which are beyond our control.

#### Scraping and waste disposal:

Dispose of the equipment and any components removed from it safely in accordance with all local and national safety and environmental requirements. Particular care must be taken with components and waste oil which have been contaminated with dangerous substances from the process. Do not incinerate fluoroelastomer seals and O-rings.

- You may authorize us to dispose of the equipment **at your expense**.

# Health and safety clearance form

## Declaration concerning safety, potential hazards and safe disposal of waste, e. g. used oil.

Safety and health of our staff, laws and regulations regarding the handling of dangerous goods, occupational health and safety regulations, safety at work laws and regulations regarding safe disposal of waste, e. g. waste oil, require that for all pumps and other products this form must be sent to our office duly completed and signed before any equipment is dispatched to our premises. **Products will not be accepted for any procedure, and handling and repair / DKD calibration will not start before we have received this declaration.**

- Fax or post a **completed copy of this form** to us in advance. The declaration must arrive before the equipment. **Enclose a second, completed copy with the product.** If the product is contaminated you must notify the carrier (**GGVE, GGVS, RID, ADR**).
- Inevitably, the repair process will be delayed considerably, if this information is missing or this procedure is not obeyed. We hope for your understanding for these measures which are beyond our control and that you will assist us in expediting the repair procedure.
- Make sure that you know all about the substances which have been in contact with the equipment and that all questions have been answered correctly and in detail.**

**1. Product (Model):** .....

**2. Serial No.:** .....

**3. List of substances in contact with the equipment or reaction products:**

**3.1 Chemical/substance name, chemical symbol:**

a) .....

b) .....

c) .....

d) .....

**3.2 Important information and precautions, e. g. danger classification:**

a) .....

b) .....

c) .....

d) .....

**4. Declaration (please mark as applicable):**

**4.1 for non dangerous goods:**

We assure for the returned product that

- neither toxic, corrosive, biologically active, explosive, radioactive nor contamination dangerous in any way has occurred.
- the product is free of dangerous substances.
- the oil or residues of pumped media have been drained.

**4.2 for dangerous goods:**

We assure for the returned product that

- all substances, toxic, corrosive, biologically active, explosive, radioactive or dangerous in any way which have been pumped or been in contact with the product are listed in 3.1, that the information is complete and that we have not withheld any information.
- the product, in accordance with regulations, has been
  - cleaned
  - decontaminated
  - sterilized

**5. Way of transport / carrier:**

.....  
Day of dispatch to VACUUBRAND:

If the paint is damaged, we wish a repaint or a replacement of parts due to optical aspects at our expense (see "Notes on return to the factory"):

- yes
- no

**We declare that the following measures - where applicable - have been taken:**

- The oil has been drained from the product.  
**Important: Dispose of according to national regulations.**
- The interior of the product has been cleaned.
- All inlet and outlet ports of the product have been sealed.
- The product has been properly packed, if necessary, please order an original packaging (costs will be charged), and marked as appropriate.
- The carrier has been informed about the hazardous nature of the goods (if applicable).

**We assure** VACUUBRAND that we accept liability for any damage caused by providing incomplete or incorrect information and that we shall indemnify VACUUBRAND from any claims as regards damages from third parties.

**We are aware** that as expressed in § 823 BGB (Public Law Code of Germany) we are directly liable for injuries or damages suffered by third parties, particularly VACUUBRAND employees occupied with handling/repairing the product.

Signature: .....

Name (print): .....

Job title (print): .....

Company's seal: .....

Date: .....

**VACUUBRAND GMBH + CO KG**

-Technology for Vacuum Systems-  
© 2008 VACUUBRAND GMBH + CO KG Printed in Germany

Alfred-Zippe-Str. 4 - 97877 Wertheim

Tel.: +49 9342 808-0 - Fax: +49 9342 808-450

E-Mail: info@vacuubrand.de - Web: www.vacuubrand.com

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Disclaimer: Our technical literature is only intended to inform our customer. The validity of general empirical values and results obtained under test conditions for specific applications depend on a number of factors beyond our control. It is therefore strictly the users' responsibility to very carefully check the validity of application to their specific requirements. No claims arising from the information provided in this literature will, consequently, be entertained.

**VACUUBRAND GMBH + CO KG**

-Vakuumtechnik im System-

© 2011 VACUUBRAND GMBH + CO KG Printed in Germany

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